

Integrated Screening in Hit Identification: Combining DEL Hit Results with FBDD and Computational Approaches

WuXi Biology


Zuyuan Shen, Zaihong Zhang, Jianyu Li, Yunjing Jia, Qi Zhang, Junkun Lei, Shuai Han, Wen Luo, Haozhen Han, Huan Liu, Jingsi Zhao, Weiren Cui, Robert Byrne, Sandra Postel, Andreas Schoop, Alex Satz, Wenji Su, Letian Kuai

Abstract

Despite the availability of various modern screening technologies and the growing experience with new modalities, poorly druggable targets still remain a challenge in the identification of hits, leads, and clinical candidates for novel targets. Often, a deeper understanding of the target and its interaction partners is required and enabling technologies like structure biology with crystal structures or cryo-EM structures, tool compounds, and specifically designed assays are key to advance projects to a stage in which medicinal chemistry can gain traction. Generating hit matter and tool compounds with one screening technology and screening with a second technology can then become quite powerful. Here, we present a case of p38 α screening in a combination of three screening technologies and structure biology using fragment screening, DEL screening, and a computational model to identify and advance hits into nanomolar small molecules.

Integration of Hit-ID Strategies & Identification of a Modality


Screening-Platforms & Chemical Space Availability

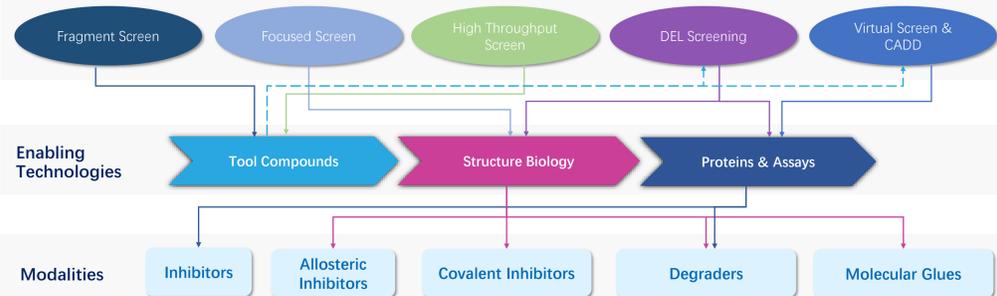

Fragment Screening & Structure Biology

Fragment Screening Workflow:

Target for p38 α (MAPK14) plays a pivotal role in initiating different disease states such as inflammatory disorders, neurodegenerative diseases, cardiovascular cases, and cancer. As a result, the identification of potent small molecule p38 α inhibitors has been actively pursued by many academic and industry groups.

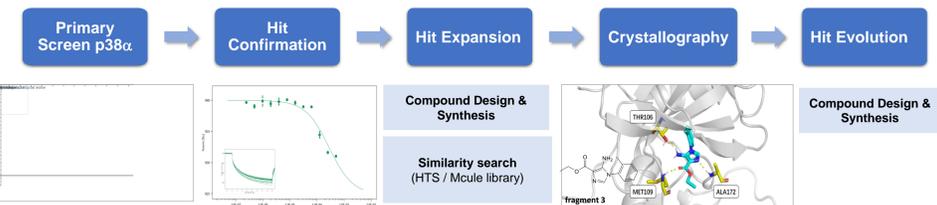


Fragment Hits Submitted to Crystallography:

Fragment	SPR K_D	LE	HAC	MW	IC_{50}	ΔT_m
Fragment 1	26.6 μ M	0.35	18	244	30.5 μ M	3.5 $^{\circ}$ C
Fragment 2	66.6 μ M	0.30	19	279	25.6 μ M	
Fragment 3	71.6 μ M	0.31	18	245		2.2 $^{\circ}$ C
Fragment 4	117 μ M	0.32	17	250	21.1 μ M	4.1 $^{\circ}$ C
Fragment 5	150 μ M	0.33	16	217		$\Delta T_m < 0.5$ $^{\circ}$ C
Fragment 6	313 μ M	0.32	15	205.31		$\Delta T_m < 0.5$ $^{\circ}$ C
Fragment 7	208 μ M	0.51	10	151	100 μ M	1.5 $^{\circ}$ C
Fragment 8	243 μ M	0.35	14	191		$\Delta T_m < 0.5$ $^{\circ}$ C
Fragment 9	365 μ M	0.43	11	250		$\Delta T_m = -1.3$ $^{\circ}$ C
Fragment 10	372 μ M	0.42	11	217		$\Delta T_m < 0.5$ $^{\circ}$ C
Fragment 11						

IC₅₀ was determined using ADP-Glo assay *ΔT_m was determined by nanoDSF*


DEL Screening

Primary hit Identified from DEL Screening

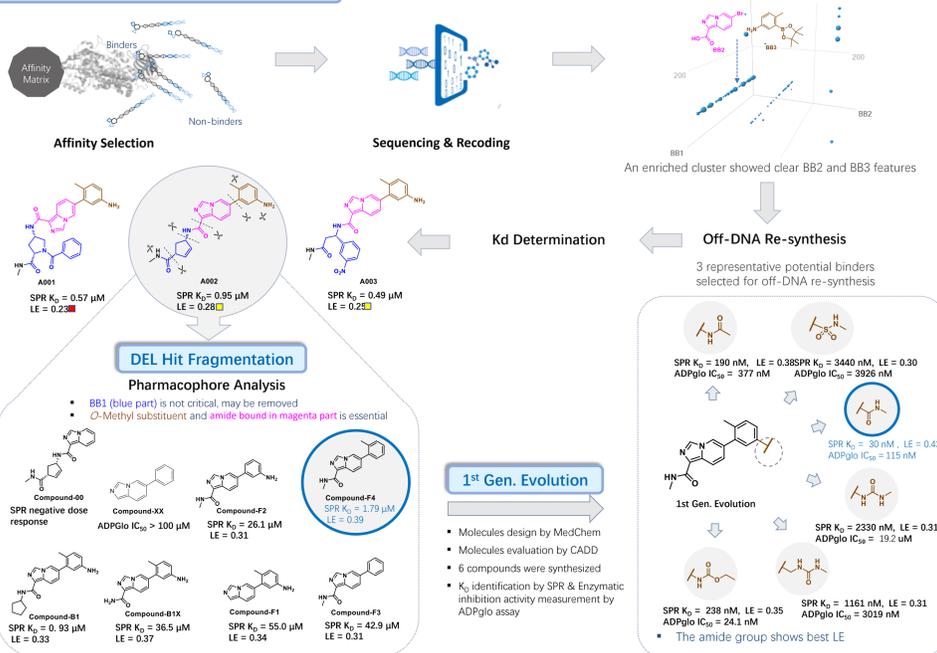

2nd Gen. Evolution


Co-crystal structure confirmation

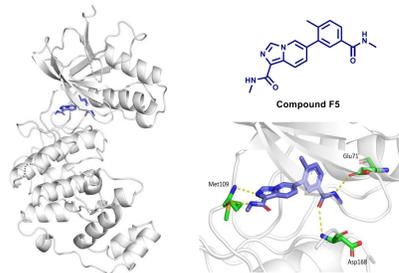

Molecule design based on co-crystal structure and MedChem knowledge

- Add aliphatic ring
- Add substituted benzene ring
- Add aromatic ring
- Add urea linked group
- Cyclization in amide
- Change the amide group to other functional groups

Optimization supported by screening of virtual molecules

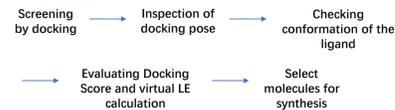

Enzymatic assay of 2nd gen. modified molecules


Compound	R ₁	p38 α IC ₅₀ (nM) ^a	LE ^b
48		81.9	0.36
16		185.0	0.33
83		12059	0.21
82		181.7	0.27
45		65.7	0.34
47		31.7	0.31
13		13.8	0.39

^a p38 α inhibition activity of compounds is tested ADP-Glo assay
^b LE value is calculated by Ki


SAR Analysis:

- Hydrophobic group is more potent than polar group for substitution
- Cyclopropyl (3-membered) substituted group shows very high ligand efficiency
- Aryl and aromatic rings are very potent


3rd Gen. Evolution

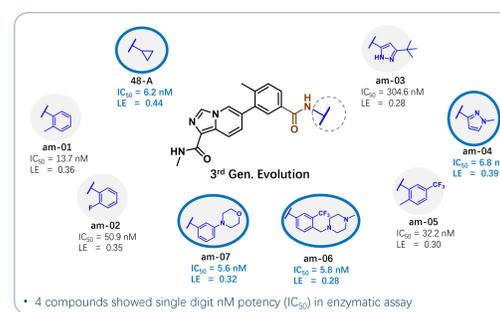

DEL Hit Fragment Evolution - Summary

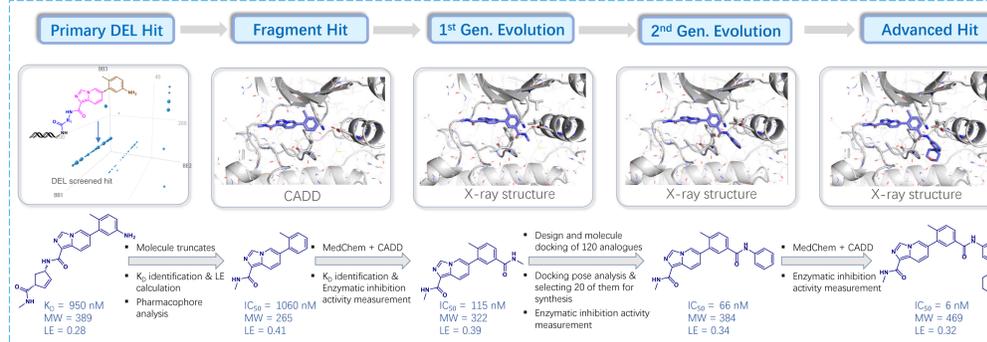

Computational Model

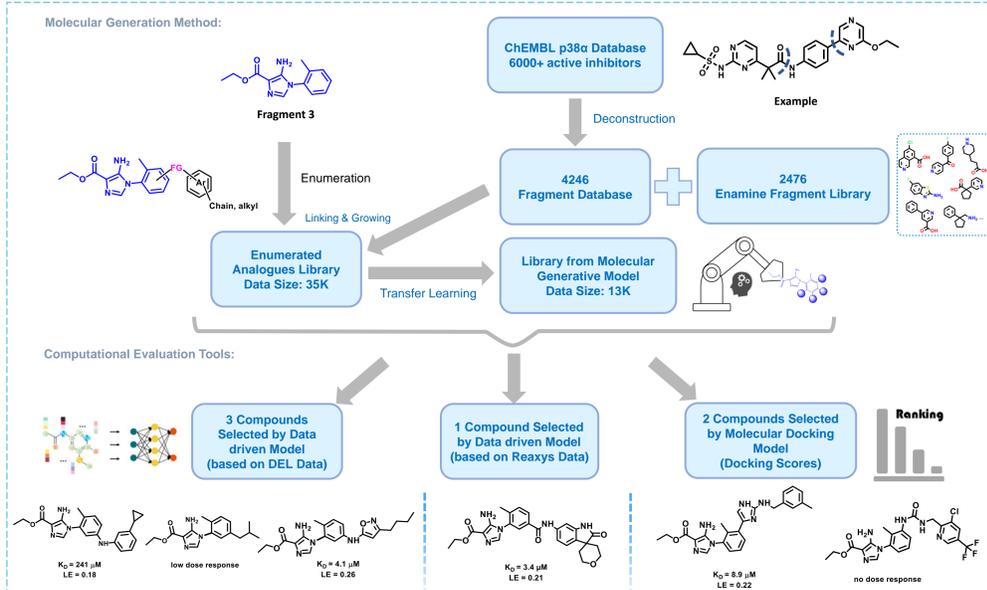

Summary

- As an integrated screening approach, we have been combining DEL-hits with fragment-based optimization methods and DEL data to improve fragment hits with computational screening approaches.
- Our DEL-screen delivered a micromolar DEL-hit that was first fragmented, and the fragment then subjected to fragment evolution to obtain nanomolar potent hits.
- Enumerating a large virtual chemical space of 35K compounds around fragment 3 derived from the fragment screen and selecting compounds for synthesis based on a DEL-model, Reaxys data or by docking resulted in micromolar potent hits.

